
### The Most Powerful Instrument

### **DESIGNED TO ADVANCE YOUR SCIENTIFIC PURSUITS**

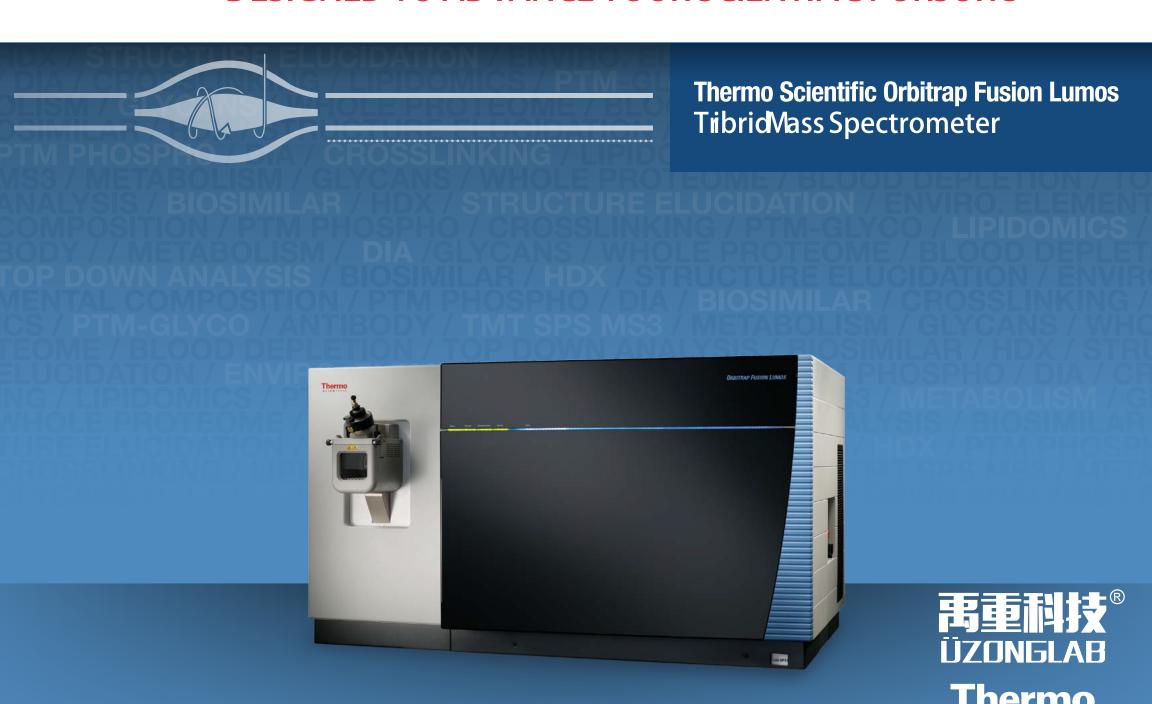

# PUSHING NEW FRONTIERS IN SCIENCE

Breakthroughs in science rely on great scientists making rapid progress answering challenging questions.

"How does lipid metabolism affect cancer?"

"How do non-cancerous cells mutate?"

"What protein modifications are indicative of a disease state?"

As a leading scientist, you need to be equipped with the latest innovations in analytical technology.

With the Thermo Scientific<sup>™</sup> Orbitrap Fusion<sup>™</sup> Lumos<sup>™</sup> Tribrid<sup>™</sup> MS, we are committed to keeping pace with your scientific pursuits through continuous innovations.



### The Most Powerful Instrument

#### Designed for High Impact Discoveries


The Thermo Scientific™ Orbitrap Fusion™ Lumos™ Tribrid™ MS delivers superior sensitivity, selectivity and versatility to enable life scientists to obtain the highest quality data. It is designed to pursue the most difficult analyses, including multiplexed quantitation of low-abundance peptides in complex matrices, characterization of positional isoforms of intact proteins, resolution of isobaric metabolites, protein structure characterization using chemical crosslinking and the deepest mining of challenging post-translational modifications.

You trusted us to make this most versatile and powerful mass spectrometer a reality. We now trust you with groundbreaking science and high impact discoveries.

## Multiplexing Targeted Analysis


#### Accurately Quantifying Hundreds of Proteins in 90 min

Multiplexed analyses using isobaric mass tags are widely utilized for high throughput quantitative comparisons of protein abundances. The TMT SPS MS³ workflow available on the Thermo Scientific™ Orbitrap Fusion™ Lumos™ Tribrid™ MS is a proven method which enables the simultaneous analysis of 10 samples together with improved quantitative accuracy. Pushing the multiplexing frontiers even further, targeted assays incorporating this technique can now be built to detect and quantify very low levels of target peptides even from undetected precursors. This new TMT SPS tMS³ workflow further benefits from the enhanced sensitivity of the Orbitrap Fusion Lumos MS, which boosts the number of quantifiable peptides present at low levels.

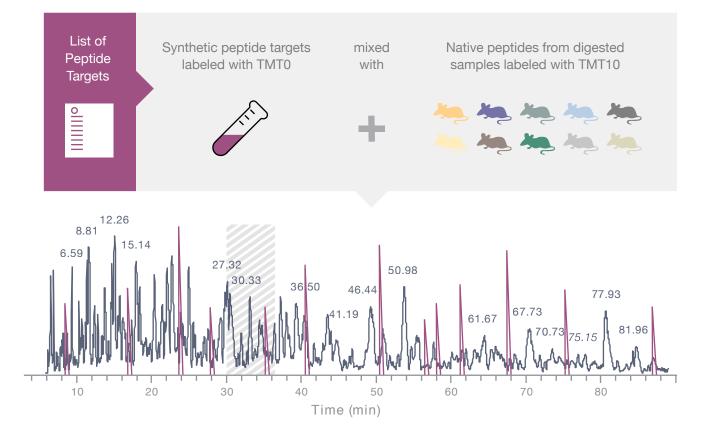

#### Base peak Extracted Ion Chromatogram (XIC). Purple peaks are XICs of TMT0-labeled peptides.

#### MS DETECTION

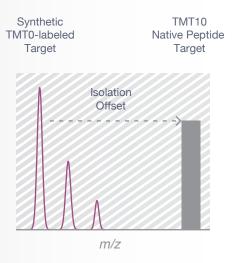


The synthetic TMT0-labeled peptide is detected in the full MS scan. The native peptide is isolated by using an offset (*m/z* difference between TMT0 and TMT10).

Everley et al. WP 663, ASMS 2016.

The Thermo Scientific™ Orbitrap Fusion™ Lumos™ Tribrid™ MS now supports targeted proteomics assays with multiplexing in two dimensions: targets and samples. These experiments make use of triggered-by-offset-in-mass peptides that allow for quantitation, even from undetected precursors. Using this method, we monitored 131 target peptides corresponding to 69 proteins across the NCI-60 cell line in biological triplicate, analyzing 180 samples in only 48 hours (16 min/sample). We found that accurate and reproducible TMT SPS tMS³ quantitation elucidated a correlation between expression of key proteins and cellular response to drug treatment.


-Steven Gygi, Professor, Harvard Medical School

#### CID MS<sup>2</sup>

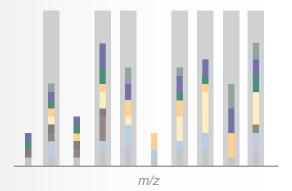


CID MS<sup>2</sup> spectrum of native peptide target contains multiplexed fragments. Gray bars indicate the predicted fragment target ions that are used for the following SPS tMS<sup>3</sup> step.

#### FTMS HCD SPS tMS<sup>3</sup>

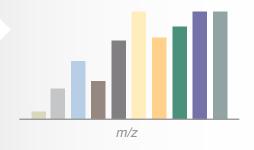

Targeted SPS MS<sup>3</sup> of the predicted MS/MS fragments generates ten reporter ions used for quantitation.

### IMPROVED QUANTITATIVE ACCURACY AT LOW LEVELS

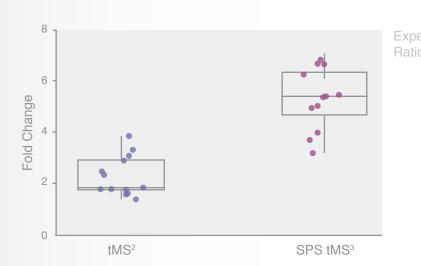



Comparison of results for the TMT tMS² and TMT SPS tMS³ quantification methods shows improved accuracy of the tMS³ method. Each point represents the fold change of a single quantification event for each of the targeted peptides present at 100 amol level in matrix. The expected ratio is 8:1.

## Advancing Peptide Quantitation

#### Large Scale Quantitation with Data Independent Acquisition


Data Independent Acquisition (DIA) is widely used today for the global identification and quantification of thousands of peptides in complex mixtures. Step-wise isolation and MS/ MS fragmentation of all ions in a defined *m/z* window cover the targeted mass range and provide ultra-high resolution MS and MS/MS data for all components in the sample, allowing accurate quantitation of peptides, and the unique opportunity for retrospective analysis of unknowns and new targets of interest in the future. The Thermo Scientific™ Orbitrap Fusion™ Lumos™ Tribrid™ MS provides speed, selectivity and enhanced sensitivity to obtain maximal performance while maintaining high reproducibility of quantitation for low abundance analytes.

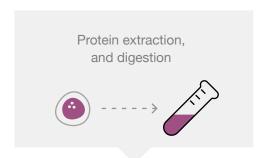

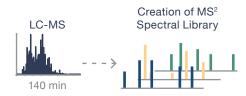


HeLa digest standard was introduced into the Orbitrap Fusion Lumos MS using RP-HPLC. Data Dependent experiments were used to generate high resolution accurate mass (HRAM) MS² spectra. The data was searched using SEQUEST™ HT in order to generate a spectral library for the following DIA data analysis.

#### DATA INDEPENDENT HRAM HCD MS<sup>2</sup>

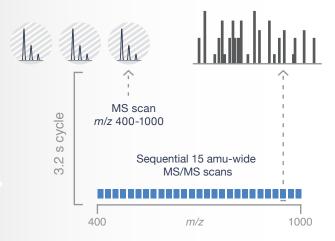

DIA acquisition was performed using a 120,000 FWHM resolution MS scan followed by 15 amu-wide MS/MS windows at 30,000 FWHM resolution with 60 ms maximum injection time.

### MATCHING DIA SPECTRUM TO THE SPECTRAL LIBRARY

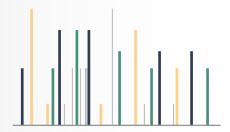


DIA MS² spectra were searched against the library built from Sequest HT search results. Fragment peak detection was performed with Biognosys Spectronaut™ software.

### QUANTITATION USING FULL MS




Peptide quantitation took into account all confidently identified isotopes from consecutive MS full scans, correcting for interference and using cross-run normalization.

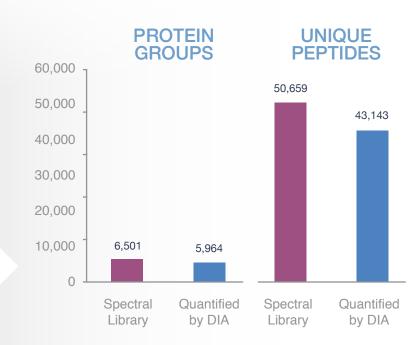

Nearly 92% of the protein groups and over 85% of the unique peptides were confidently matched to the spectral library and quantified with <6.9% CV when using 500 ng of HeLa digest.

### Confident Low Attomole Limit Peptide Quantitation Using Parallel Reaction Monitoring

Parallel Reaction Monitoring (PRM) is uniquely designed for quantifying hundreds of targeted proteins in complex matrices. Using this approach, precursor ions are isolated and selectively fragmented, with the resulting product ions analyzed in the Orbitrap. This approach benefits from the brighter ion source and Advanced Quadrupole Technology of the Thermo Scientific™ Orbitrap Fusion™ Lumos™ Tribrid™ MS, routinely achieving attomole-level limits of quantification (LOQ) in matrix.

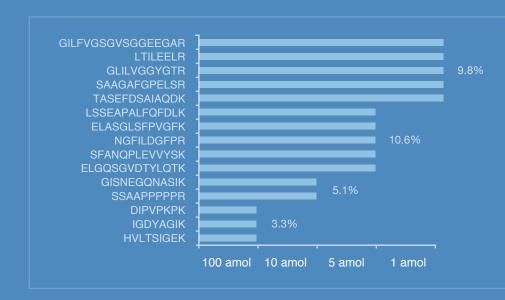

Fifteen (15) PRTC peptides were spiked into 200 ng of HeLa digest and analyzed by LC-MS (30 min run). The Orbitrap Fusion Lumos MS provides accurate quantitation of all 15 PRTC peptides with some down to 1 attomole levels. Average CV% for each LOQ level is shown.

## New Horizons in Intact Protein Analysis


#### Highly Selective Analysis of Protein Isoforms

Top-down mass spectrometry is commonly utilized to characterize intact proteins and their modifications. The Advanced Vacuum Technology, unique to the Thermo Scientific™ Orbitrap Fusion™ Lumos™ Tribrid™ MS, provides conditions optimized for improved performance in intact protein analysis. The high selectivity of Advanced Quadrupole Technology allows for isolation of precursors and detection of fragments with very high resolving power in the Orbitrap analyzer. Combined, the new system most efficiently delivers the highest quality data for the characterization of protein isoforms and their post-translational modifications.

"The Lumos proved capable of high selectivity precursor selection in MS but retained exceptionally high sequence coverage for MS<sup>2</sup>. This enabled the decoding of the most complicated core histone H3 to detect trivalent proteoforms uniquely detected in a cellular model of the B-cell cancer, multiple myeloma."

-Neil Kelleher, Professor, Northwestern University


### SPECTRUM, 18+

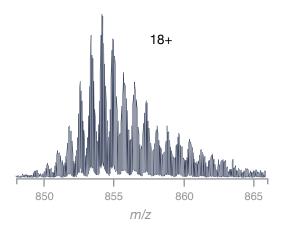

Advanced Vacuum Technology increases ion transmission, improving the quality of intact protein spectra acquired in the high resolution accurate mass Orbitrap analyzer.

Zheng, Y et al. Mol Cell Proteomics 2016, 15,776-90.

## HIGH RESOLUTION ISOLATION OF METHYLATED FORMS OF HISTONE H3

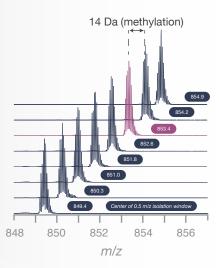

With improved ion transmission provided by the Advanced Quadrupole Technology, it is now possible to efficiently enrich for individual isobaric protein forms for a subsequent top-down analysis.

## FRAGMENTATION MAPS OF HISTONE H3 FROM ETD HD™ TOP-DOWN EXPERIMENTS


ETD HD identification of isomeric forms of histone H3, differing in the site of trimethylation (K9 vs. K27) from the precursors at m/z 853.4.

ETD HD enhances the dynamic range of ETD spectra by increasing the precursor ion storage capacity. The higher efficiency of ETD HD experiments provides greater sequence coverage at faster acquisition rates.

#### Comprehensive Characterization of Intact Monoclonal Antibodies

Recent advances in protein and antibody-based therapeutics have led to a demand for mass spectrometers capable of comprehensive characterization of antibody heterogeneity in addition to providing accurate mass measurements. The Thermo Scientific™ Orbitrap Fusion™ Lumos™ Tribrid™ MS allows high accuracy mass analysis of the intact monoclonal antibody with isotopic resolution of the heavy and light chains. The combination of the various fragmentation methods available on the Orbitrap Fusion Lumos MS and the improved efficiency provided by ETD HD enables users to obtain high sequence coverage for the light and heavy chains.

#### **INTACT mAb GLYCOFORMS**

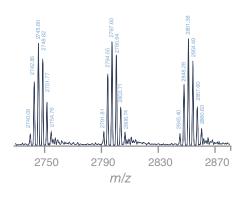



148541.91	7.0	G1F/G2F	38.6165
148380.82	0.0	G1F/G1F	84.1024
148219.21	3.6	G0F/G1F	100.0000
148058.29	11.8	G0F/G0F	68.152
147909.08	9.2	G0/G0F	10.2611

Average Matched Matched Relative
Mass Delta Glycoforms Abundance
Mass (ppm)

Raw spectrum showing baseline resolution of major intact mAb glycoforms. Deconvolution results are shown in the table above.

### ISOTOPIC RESOLUTION OF mAb LIGHT AND HEAVY CHAINS

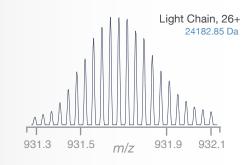

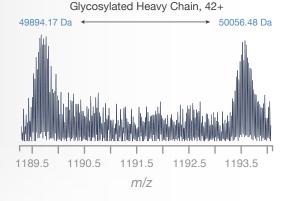

Isotopic resolution of light chain (top panel), and heavy chain (bottom panel), showing the detection of glycosylated forms.

### TOP-DOWN SEQUENCE COVERAGE FOR mAb LIGHT AND HEAVY CHAINS

```
N D V(L M\T\Q\T\P\L\S\L\P\V\S\L\G\D\Q A S I S C\R\S 25
26\S\Q\Y\I\V H\S\N\G\N\T\Y L\E\W\Y\L Q\K P\G\Q\S P\K 50
51\L\L\I\Y\K\V\S\N\R\F\S\G\V P\D\R\F\S\G\S\G\G\S
51\L\L\I\Y\K\V\S\N\R\F\S\G\V P\D\R\F\S\G\S\G\G\S\G\T\D 75
76\F\T\L\K\I\S\R\V\E\A\E\D\L\G\V\Y\Y\C\F\Q G\S\H\V P 100
101\L\T\F\G\A\G\T\K\L\E\I\K\R\A\D\A\A\P\T\V\S\I\F\P\P 125
126 S\S\E\Q\L\T\S\G\G\A\S\V\V\C\F\L\N\R\F\Y\P\K\D\I\N 150
151\V\K\W\K\I\D\G\S\E\R\Q\N\G\V\L\N\S\W\T\D\Q\D\S\K\D 175
176\S\T\Y\S\M\S\S\T\L\T\L\T\K\D\E\V\E\R\H\N\S\Y\T\C\E 200
201\A\T\H\K\T\S\T\S\P\I\V\K\S\F\N\R N E\C
```

N Q V Q L|K|E|S|G|P|G|L|V|A|P|S|Q|S|L|S|I|T|C|T|V|S 25 26 G F)SLLLGYG V NW VRQPP GQ GLEW LMG 50 51 I W G D G S T D Y N S A L K S R I S I T K D N S K 75 76 S Q V F L K M N S L Q T D D T A K Y Y C T R A P Y 100 101 GKQYFAYWGQGTLVTVSAAKTTPP S 125 126 VIYIP LIAIPIGISIAIAIOITIDIS M VIT L G C L VIKIG Y 150 151 FPEPVTVTWNSGSLSSGVHTFPAVL 175 176 O)S D L Y T L S S S V T V P S S T W P S E T V T C 200 201 N V A H P A S S T K V D K K I V P R D C G C K P C 225 226 I CTVP E VSS V F I F P P K P K D V L T I T L 250 251 T P[K V T C]V V V D I S K D]D P] E V Q[F[S|W[F V]D 275 276 DV EVHTAHTQPREEQFNSTFRSVSE 300 301 LLP I MHQDWLNGKEFK CR VN S A A F PA 325 326 PLILE KITLISKIT K GRPKA PQ V YLTLI P PPK 350 351 ELQMAKDKVSL T CM ILTDEFFPEDITTVE 375 376 WQWNGQPAENYKNTQPIMDTDGSYF 400 401 VYSKLINVQKSNWEAGNTFTCSVLLHE 425 426 GLL HINHHITLE KISLLS H S P G C

Combination of ETD HD, CID and HCD fragmentation modes provided 91% sequence coverage for the light chain (top panel) and 62% sequence coverage for the heavy chain (bottom panel).

### Revolutionizing Glycoproteomics

#### Simplified Glycopeptide Analysis

Analysis of post-translational modifications (PTMs) is often central to biological research because of their roles in cellular function and disease states. Some PTMs, such as glycosylation, are particularly challenging to analyze because they require information to both accurately determine the composition of the modification as well as correctly identify the modification site. The multiple fragmentation capabilities (CID, HCD, ETD and EThcD) on the Thermo Scientific™ Orbitrap Fusion™ Lumos™ Tribrid™ MS are essential for high-throughput, comprehensive glycopeptides analysis.

The high dynamic range of ETD and EThcD HD yields high quality MS/MS spectra of glycopeptides backbone fragmentation, resulting in higher sequence coverage and accurate localization of the modification site. Accompanying HCD spectra provide complementary information about glycan composition. Combination of fragmentation techniques, in addition to intelligent instrument control workflows for glycopeptide analysis unique to the Orbitrap Fusion Lumos MS, enables deeper mining of glycosylation modifications in a variety of biological samples.

## ORBITRAP FUSION LUMOS CAN PRODUCE:


Large scale LC-MS intact glycopeptide identification

Intelligent acquisition strategies for complete characterization of intact glycopeptide structure using HCDproduct ion dependent-EThcD/CID

Increased peptide sequence coverage and confidence in glycosylation site localization with EThcD

Quantitation of intact glycopeptides using isobaric tags via SPS MS<sup>3</sup>


#### EThcD MS/MS spectrum of O-linked glycopeptide

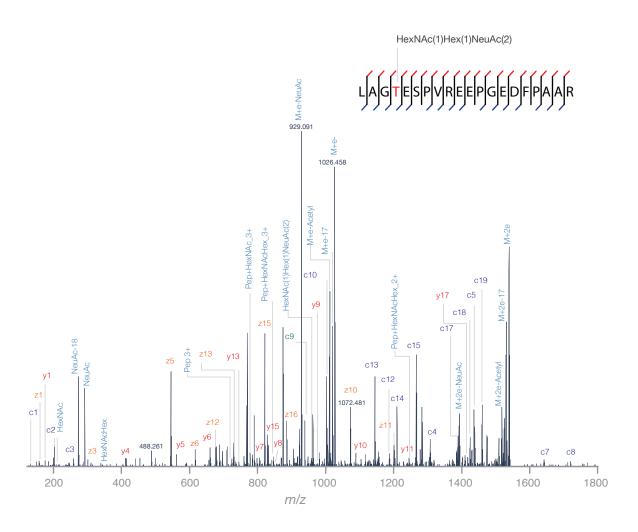

## Comprehensive Lipid Profiling

#### Enhancing Productivity with Fast Polarity Switching


For specific lipid sub-classes such as phosphatidylcholine (PC), both positive and negative HCD MS/MS data are required for full characterization of the individual molecular species. The Thermo Scientific™ Orbitrap Fusion™ Lumos™ Tribrid™ MS allows fast polarity switching and intelligent acquisition of HCD and CID MS<sup>n</sup>. This provides complementary fragmentation information for identifying PC molecular species and elucidating co-eluting triacylglycerol (TG) isomers within a single LC-MS run, yielding higher throughput, increased sensitivity and more confident analysis of lipid molecular species.

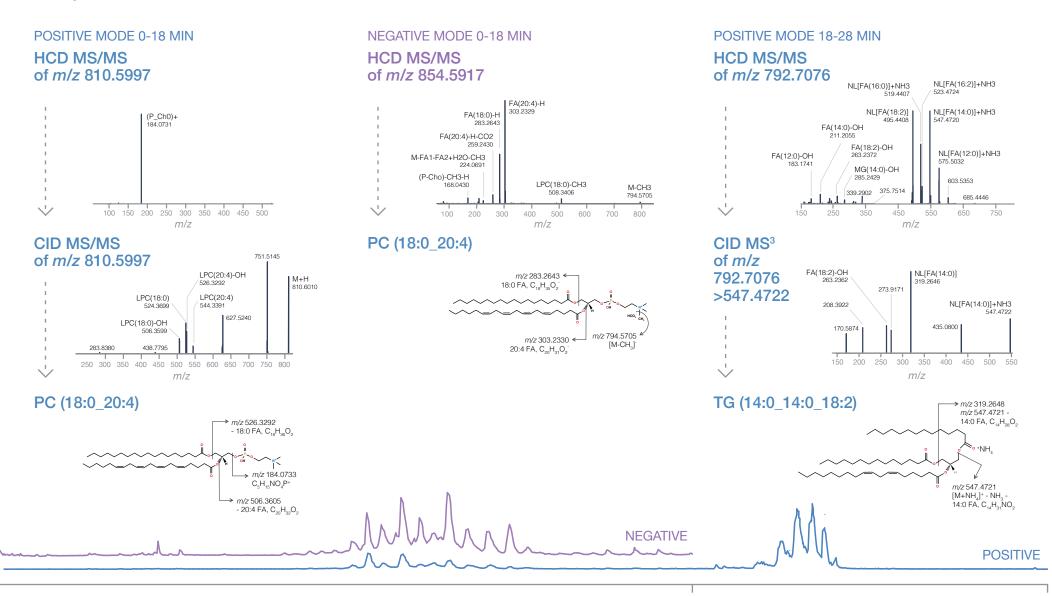


## Small Molecule Analysis


#### Ultimate Confidence in Identification and Quantitation

The high sensitivity and high resolution of the Thermo Scientific™ Orbitrap Fusion™ Lumos™ Tribrid™ MS makes it a powerful instrument for the identification and quantitation of small molecules. Due to the instrument's very high resolution capabilities, fine isotopic structure can be observed, enabling the determination of highly accurate molecular formulae. This direct measurement removes the ambiguity of pattern matching estimations and is critical in cases where monoisotopic elements like fluorine or phosphorous may be present in the compound, as shown in the example of norfloxacin below. Furthermore, the enhanced sensitivity of the Orbitrap Fusion Lumos MS enables far lower levels of quantitation.

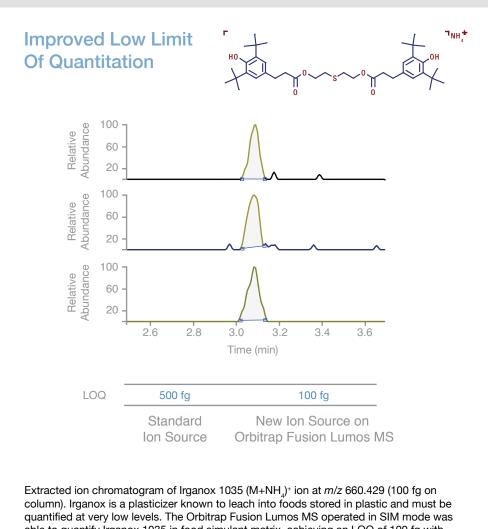

able to quantify Irganox 1035 in food simulant matrix, achieving an LOQ of 100 fg with linear dynamic range of 5 orders and <10% CV for all levels. This LOQ is 5x lower than achieved earlier on an Orbitrap instrument equipped with a standard ion source.

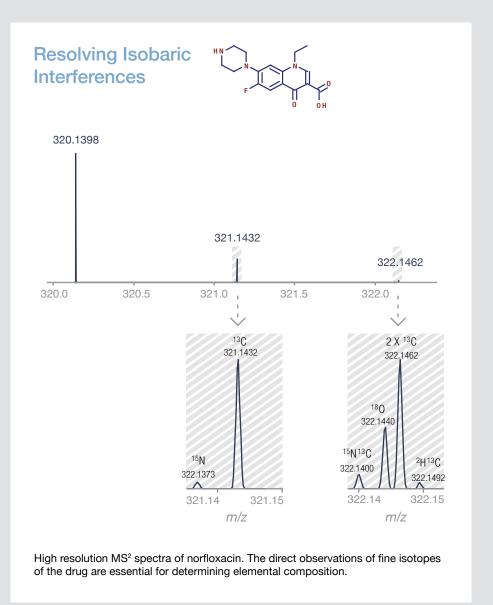

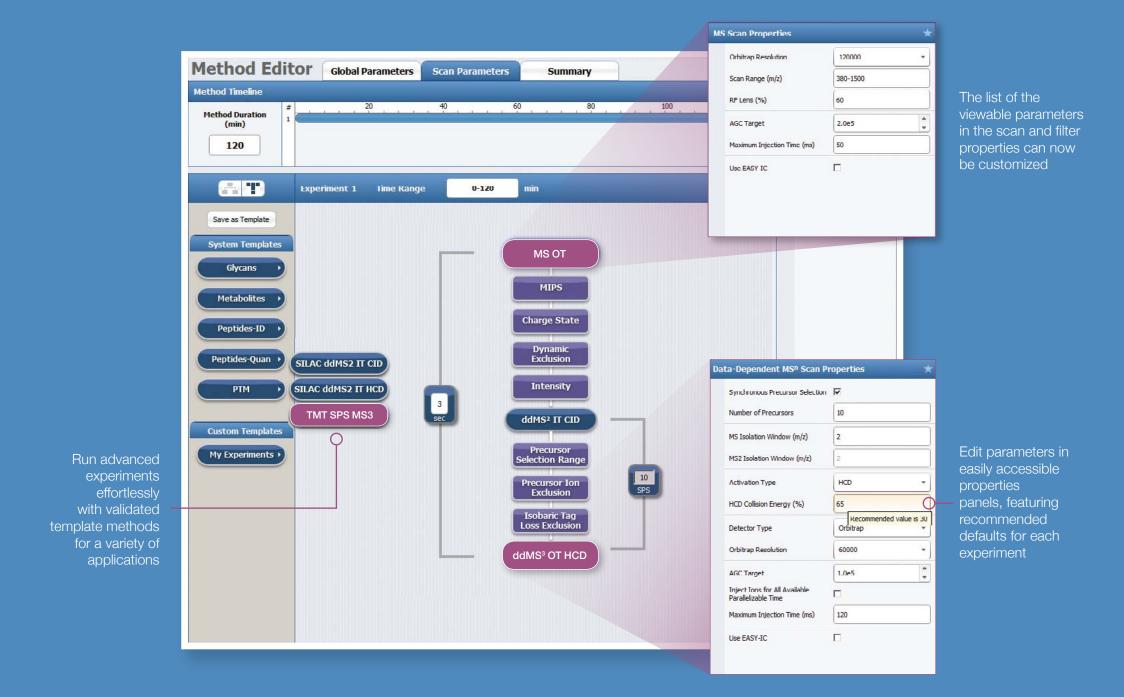

### Focus on Your Science, Not on Instrument Set Up

The highly intuitive method editor features a user-friendly interface that includes optimized pre-designed templates for a wide range of applications.



## Building on Revolutionary Tribid Architecture

The Thermo Scientific™ Orbitrap Fusion™ Lumos™ Tribrid™ mass spectrometer further amplifies the power and versatility of the innovative Tribrid design, first pioneered on the Orbitrap Fusion MS. The Orbitrap Fusion Lumos mass spectrometer incorporates the latest technologies and groundbreaking innovations in ion transmission, dissociation and detection. The combination of these improvements makes it the most sensitive, most selective and most versatile mass spectrometer to date.

With the Orbitrap Fusion Lumos MS, we are true partners, committed to innovations to advance your scientific pursuits.

### ADVANCED ACTIVE ION BEAM GUIDE


Prevents neutrals and high velocity clusters from entering mass resolving quadrupole

#### EASY ETD SOURCE

Based on Townsend discharge; reliable and easy to use


## ELECTRODYNAMIC ION FUNNEL

Focuses ions after High Capacity Transfer Tube; broad tuning curves

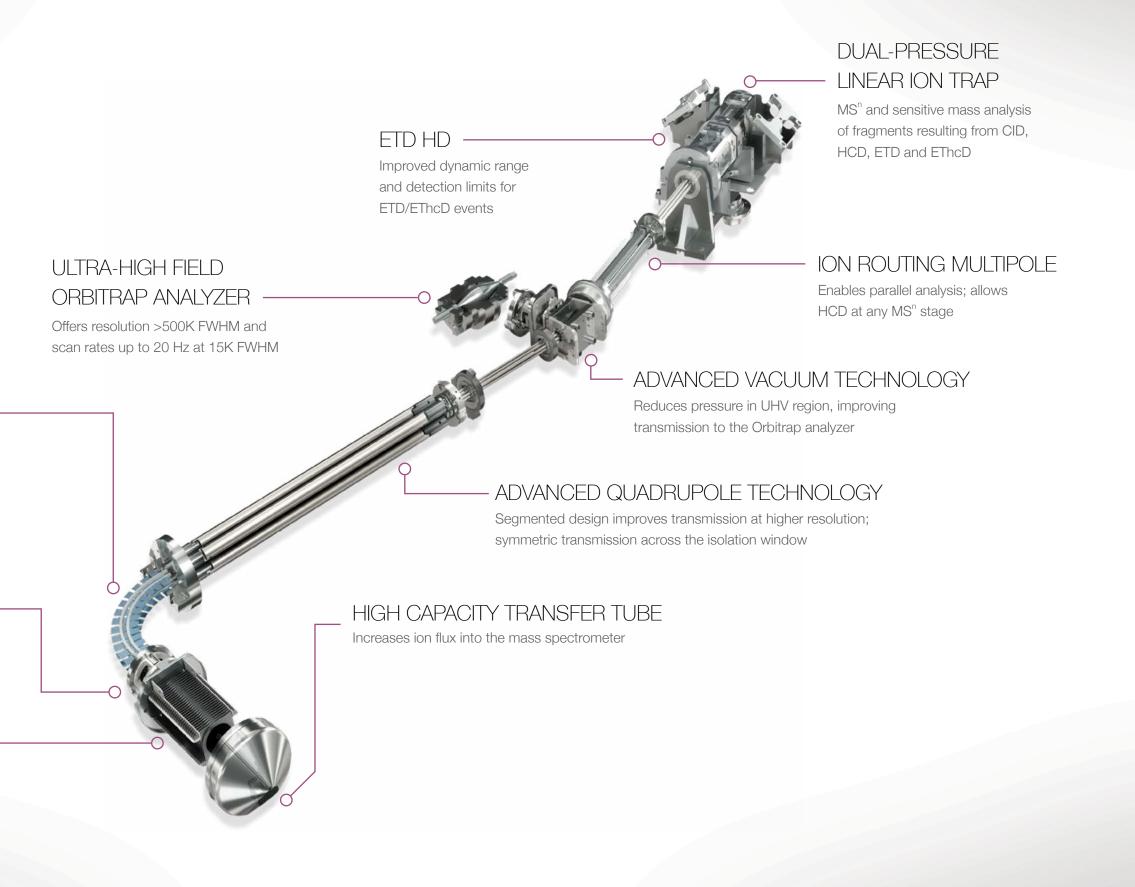

### Pushing New Frontiers in Science

The Industry's Leading Portfolio of Mass Spectrometry Solutions

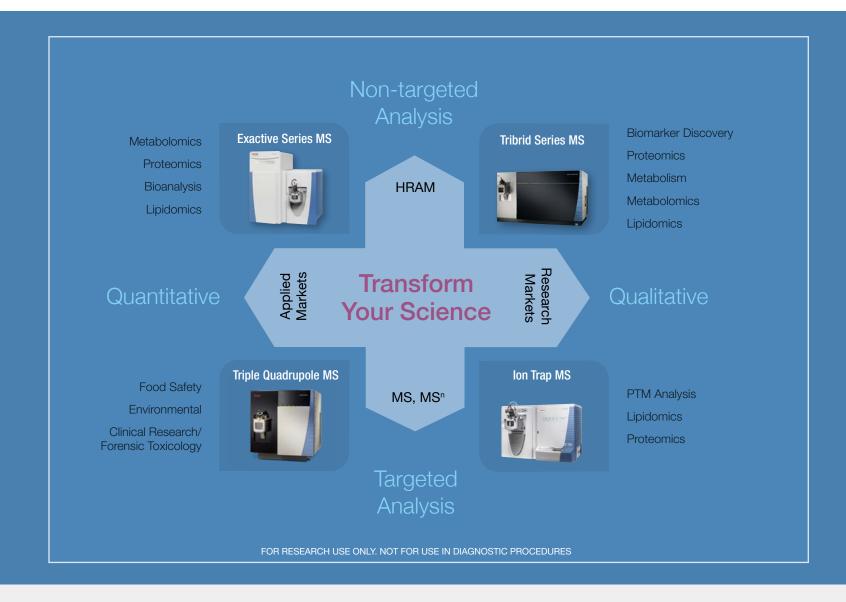

©2016 Thermo Fisher Scientific Inc. All rights reserved. SEQUEST HT is a registered trademark of the University of Washington. Biognosys Spectronaut is a registered trademark of BiognoSYS AG. All other trademarks are the property of Thermo Fisher Scientific and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

Africa Australia Austria Belgium Canada

+43 1 333 50 34 0 +61 3 9757 4300 +43 810 282 206 +32 53 73 42 41 +1 800 530 8447 800 810 5118

400 650 5118

(free call domestic)

+45 70 23 62 60 Denmark Europe-Other +43 1 333 50 34 0 +358 10 3292 200 Finland +33 1 60 92 48 00 France Germany +49 6103 408 1014 +91 22 6742 9494 India

+39 02 950 591 Italy Japan +81 45 453 9100 +82 2 3420 8600 Korea Latin America +1 561 688 8700 Middle East +43 1 333 50 34 0 Netherlands +31 76 579 55 55

New Zealand +64 9 980 6700 Norway +46 8 556 468 00 +43 1 333 50 34 0 Russia/CIS +65 6289 1190 Singapore Spain +34 914 845 965 Sweden +46 8 556 468 00

Switzerland +41 61 716 77 00 +44 1442 233555 USA +1 800 532 4752 BR64392-EN 0716M

